

# Empty bottle – Z13 & Z14

When a chemical bottle is empty please follow this procedure:

1. Rinse the inside:
  - Fill half of the bottle with DI water, close it and shake it. Empty the bottle in the sink for gloves rinsing. **Repeat the operation 3 times.** Close the bottle.
  
2. Rinse the outside:
  - Wash the outside of the bottle with a moist tissue. Then rinse the tissue with DI water. **Repeat the operation 3 times.** Or rinse the bottle 1 min in the sink for gloves rinsing.
  
3. Dry the outside:
  - Use papers to dry the outside of the bottle.
  
4. **Write on the bottle “Rinsed” + the date + your CMi username.**
  
5. Put the bottle in the grey box “FOR EMPTY AND RINSED BOTTLES” (Available in the room where you are!).

*Don't forget that people who will handle the bottle outside of the clean room will do it **without any glove** so the bottle must be very clean and dry!*