


III. Creating New Map Patterns (Quick mode)


1. Open “RSMAPPER”

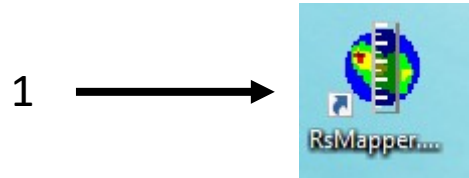

2. Click the “Wafer Map” tab



3. Select “Edit> Map Pattern” by clicking on the drop-down menu

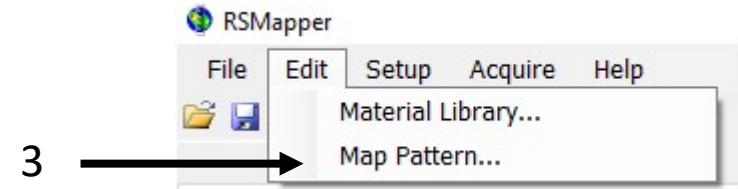


NB: The pattern that you will define in this menu is without dimension !

Your pattern will be adapted to the dimension of your sample defined in the menu


“Edit Recipe > Wafer Map” (see Editing Film Recipe)

- Set the shape of you sample ①
- Select the type of pattern ②
- Set the number of point on the ring ③
- Set the number of points on the first ring of the quadrant ④
- Set the additional point on the next ring same quadrant ⑤
- The diameter and the number of points are informative ⑥

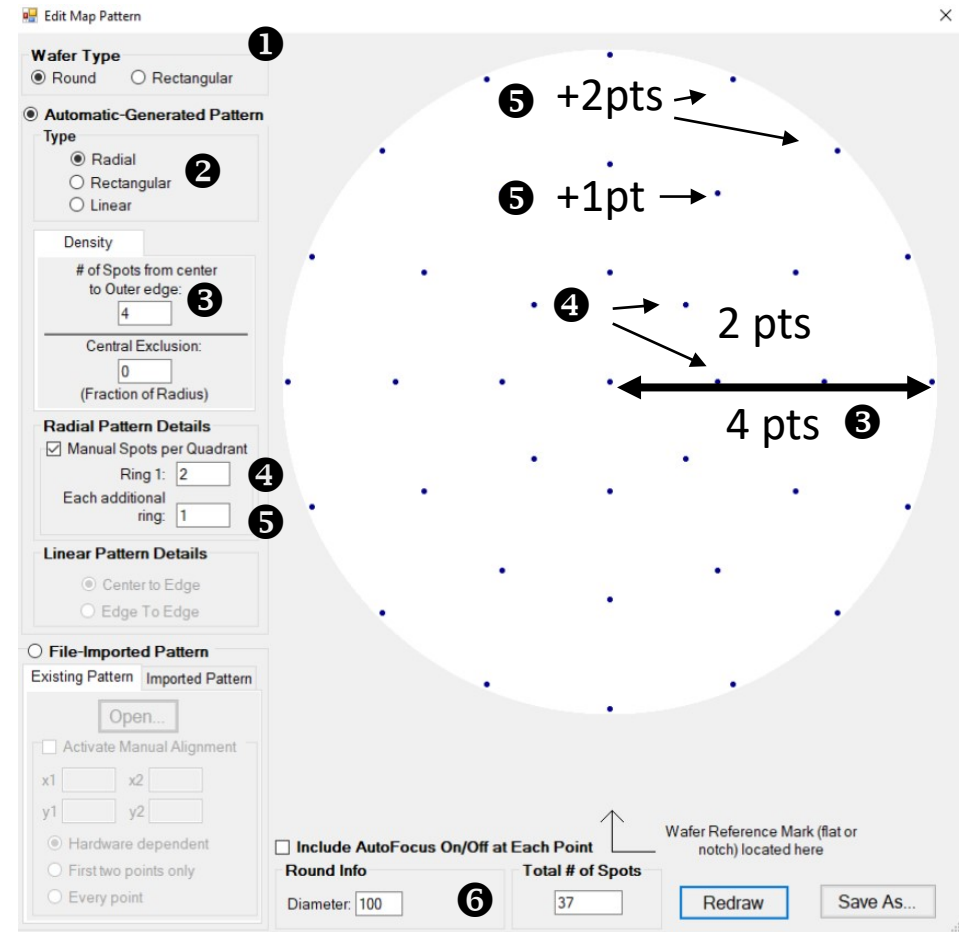


To visualize the pattern if you modify any number in the “Density” menu, press “Redraw” ↗

To save your pattern, press “Save As”

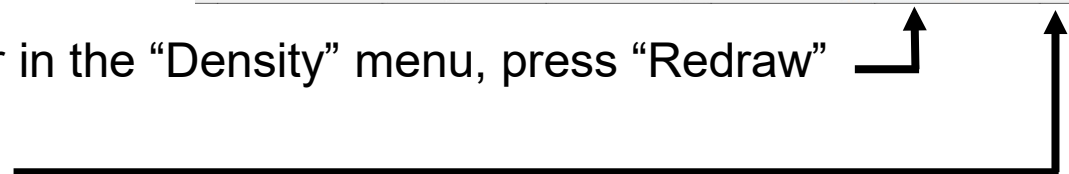

- Set the shape of you sample ❶
- Select the type of pattern ❷
- Set the number of points from the center to the outer edge ❸
- The Width & Height and the number of points are informative you can not modify them ❹

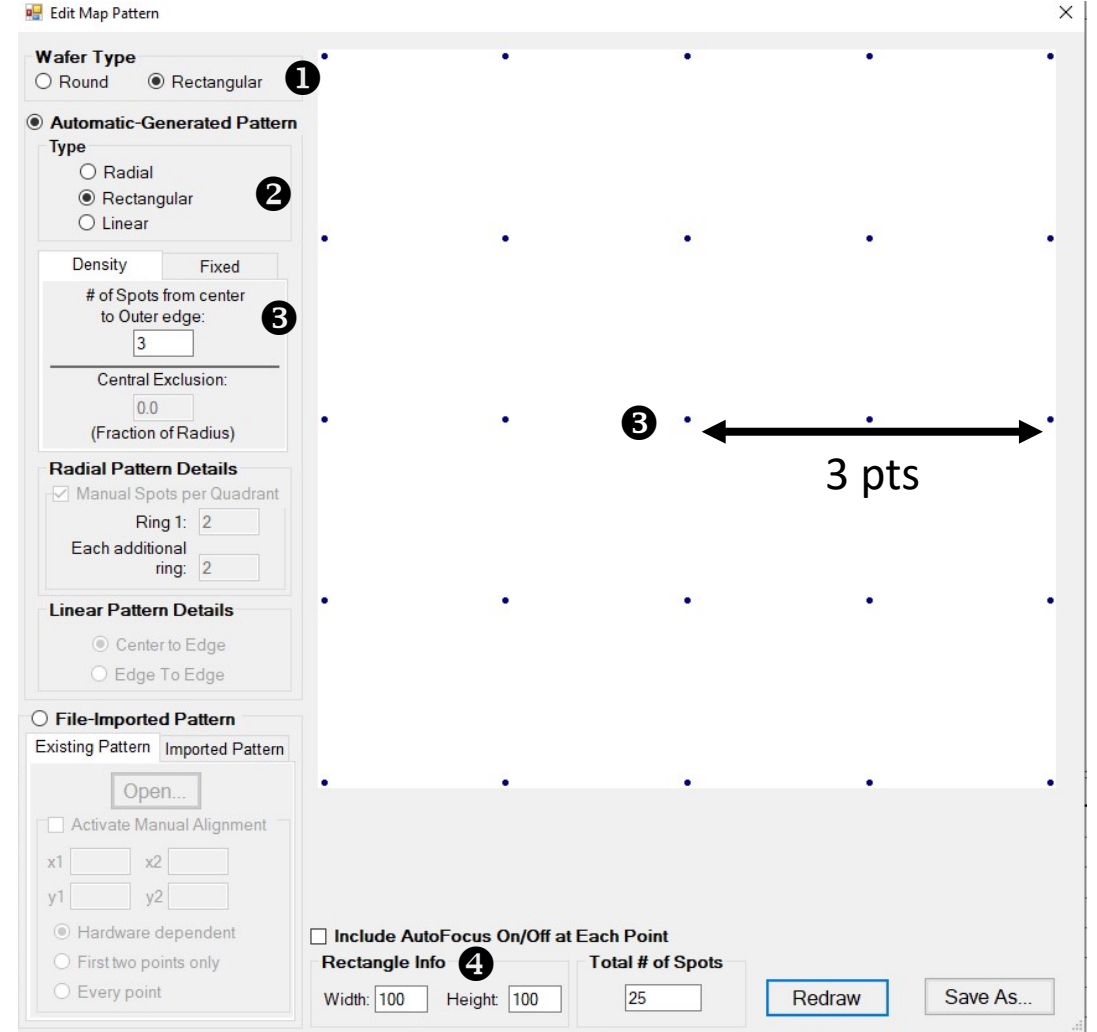

To visualize the pattern if you modify any number in the "Density" menu, press "Redraw" ↗

To save your pattern, press "Save As" ↘

