

II. Editing Film Recipe (Quick mode)


1. Open “RSMAPPER”

1 →




2. Click the “Wafer Map” tab

2 ↓

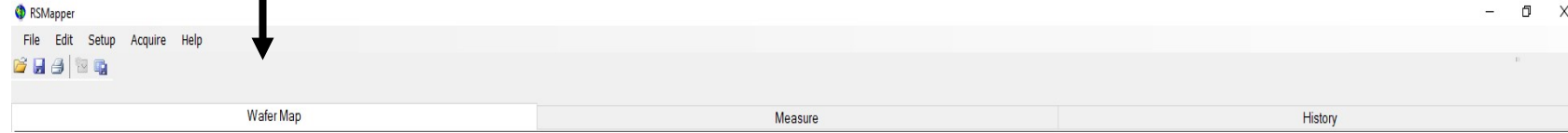


3. Select the recipe “new recipe” in the folder “CMI” by clicking on the drop-down menu

3 ↙ ↘

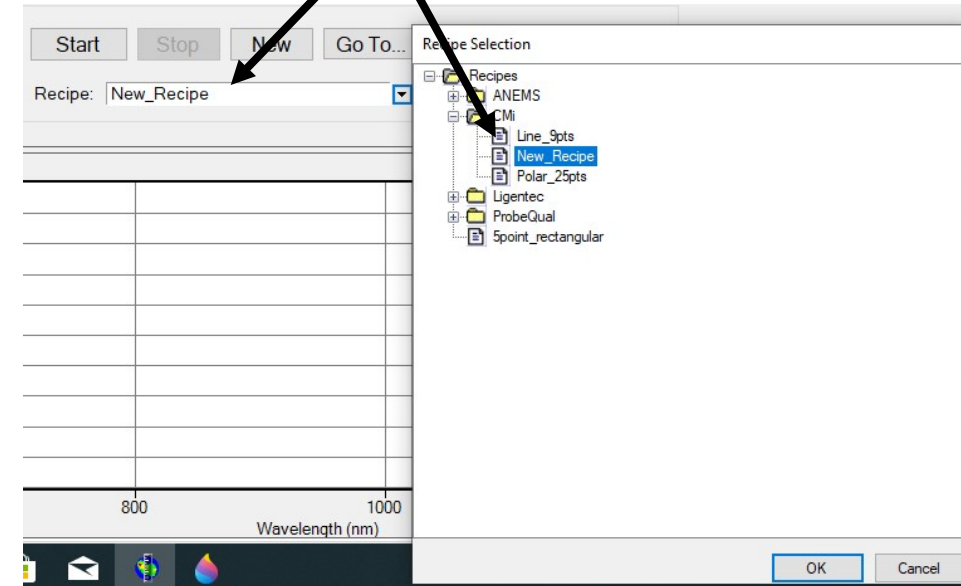

4. Click on “Edit” button

4 ↓

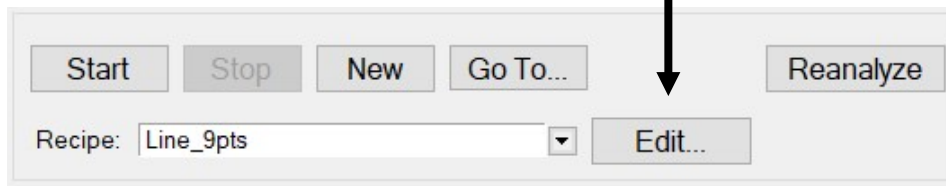

Click the “Wafer Map” tab ❶


- Define the shape of your samples ❷
Circular (Flat or notch) or rectangular
- Set the wafer diameter or the size of your sample ❸
- Select a pre-defined coordinate system with the drop-down menu ❹
- Select a pre-defined number of points with the drop-down menu ❺
- Set the edge exclusion ❻
- Set the rotation angle ❼

The screenshot shows the 'Edit Recipe' dialog box with the 'Wafer Map' tab selected. The 'Measurement Configuration' section is highlighted, and several fields are annotated with circled numbers 1 through 6. The 'Indexing' section is also annotated with a circled number 2. The 'Wafer Settings - All Sites' section is visible on the right. The 'OK', 'Cancel', and 'Apply' buttons are at the bottom right.

You can visualize the pattern once you click on “OK” or “Apply” (see “Making Measurement”)

Click the “Rs Acquisition Settings” tab ①

- Check that you are in the “Auto range” mode ②
- No need to modify the other parameters

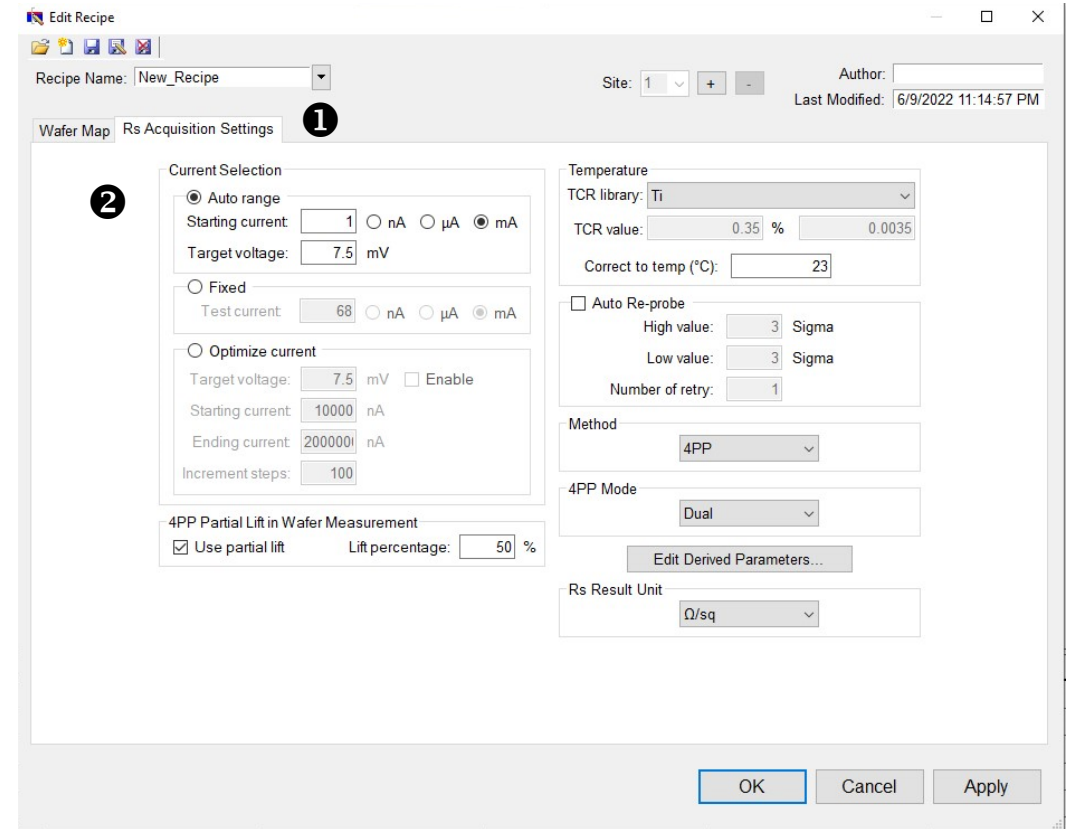

5. Save your recipe with a new name

