



I. Making Measurement (Quick mode)

1. Open “RSMAPPER”

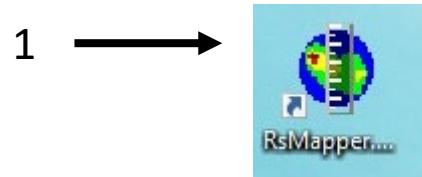


1 →

2. Click the “Measure” tab

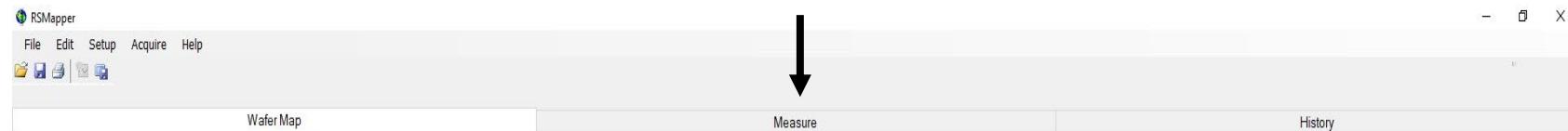


2 ↓

3. Click the “Go To” button

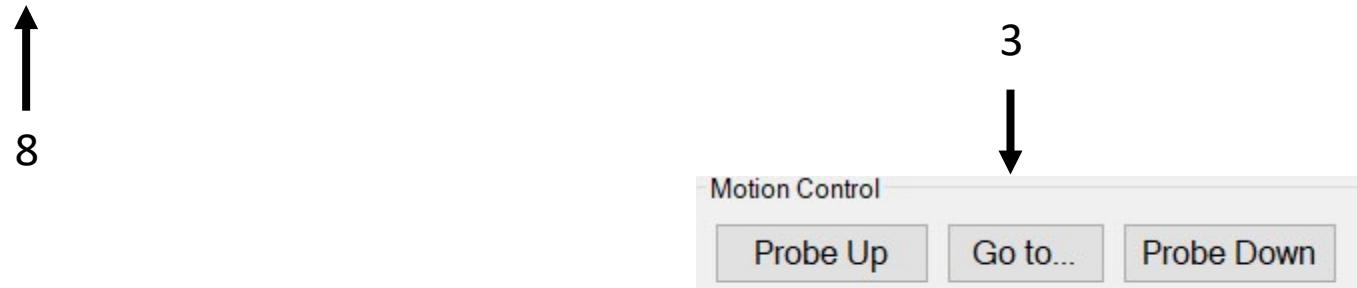

3 ↓

4. Click “Go To Unload”

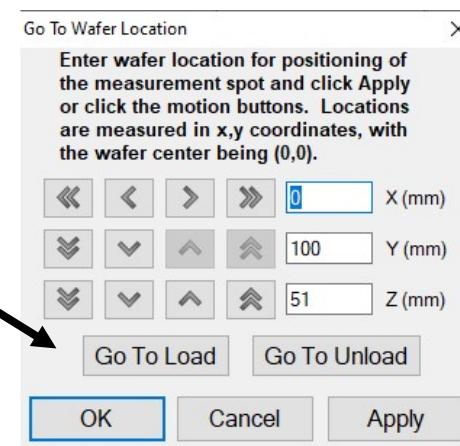

6

4

5. Load your sample/wafer

6. Click “Go To Load” and “OK”

7. Click the “Go To Z Measure”



8. Click the “Wafer Map” tab



7 ↑


9. Select your recipe by clicking on the drop-down menu

10. Click “Start”

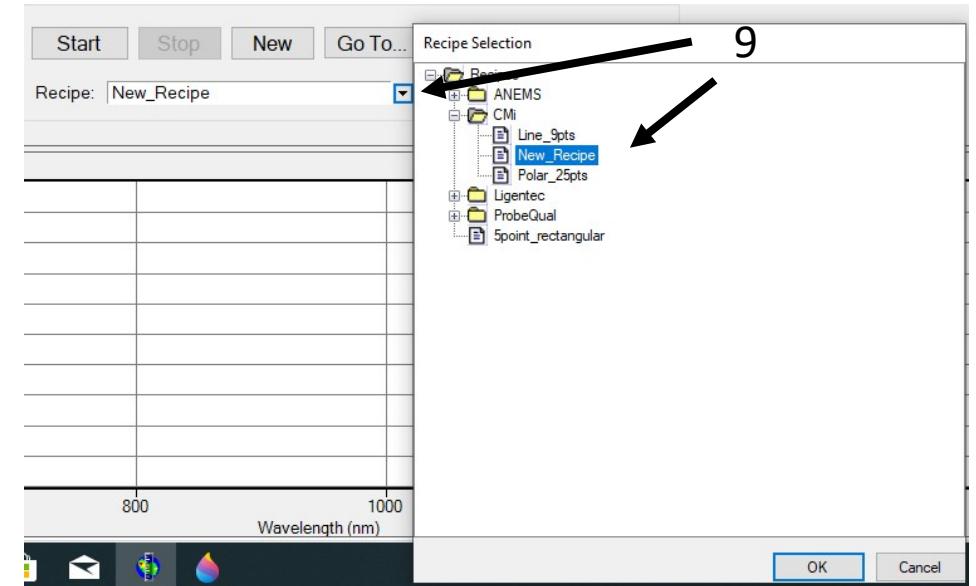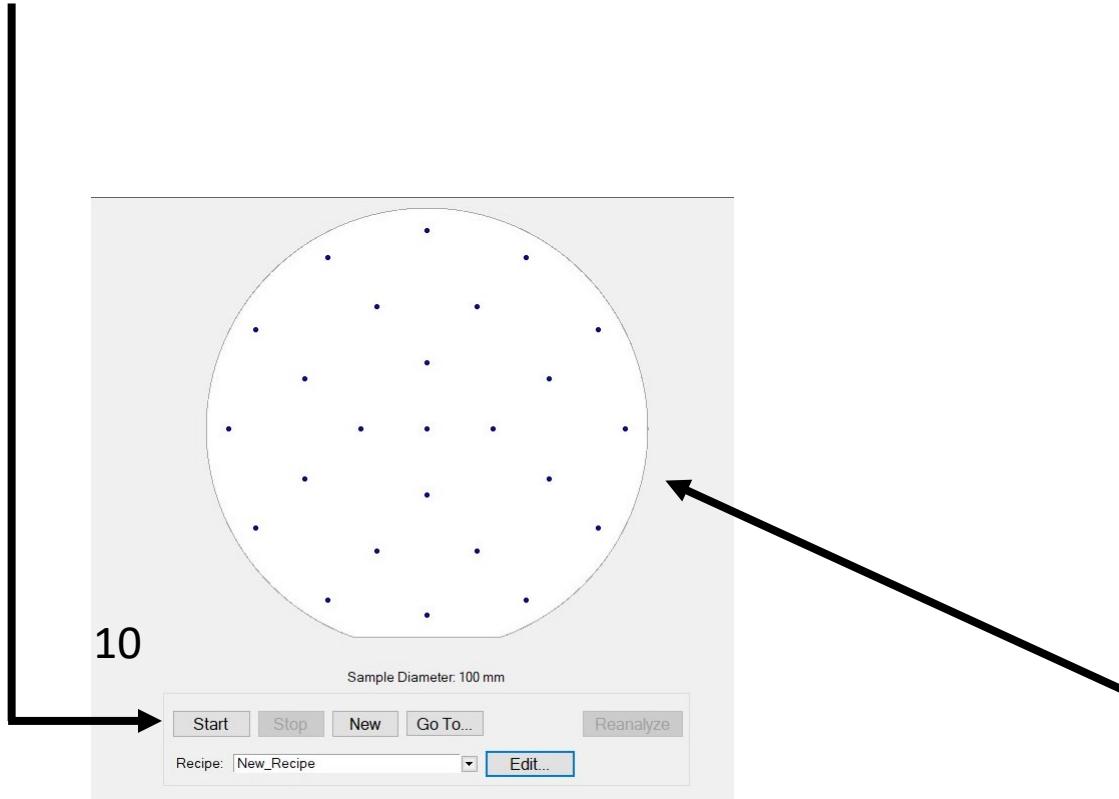

Visualization of the pattern generated by the Recipe Editor


Visualization of the measurements on the “Wafer Map” tab

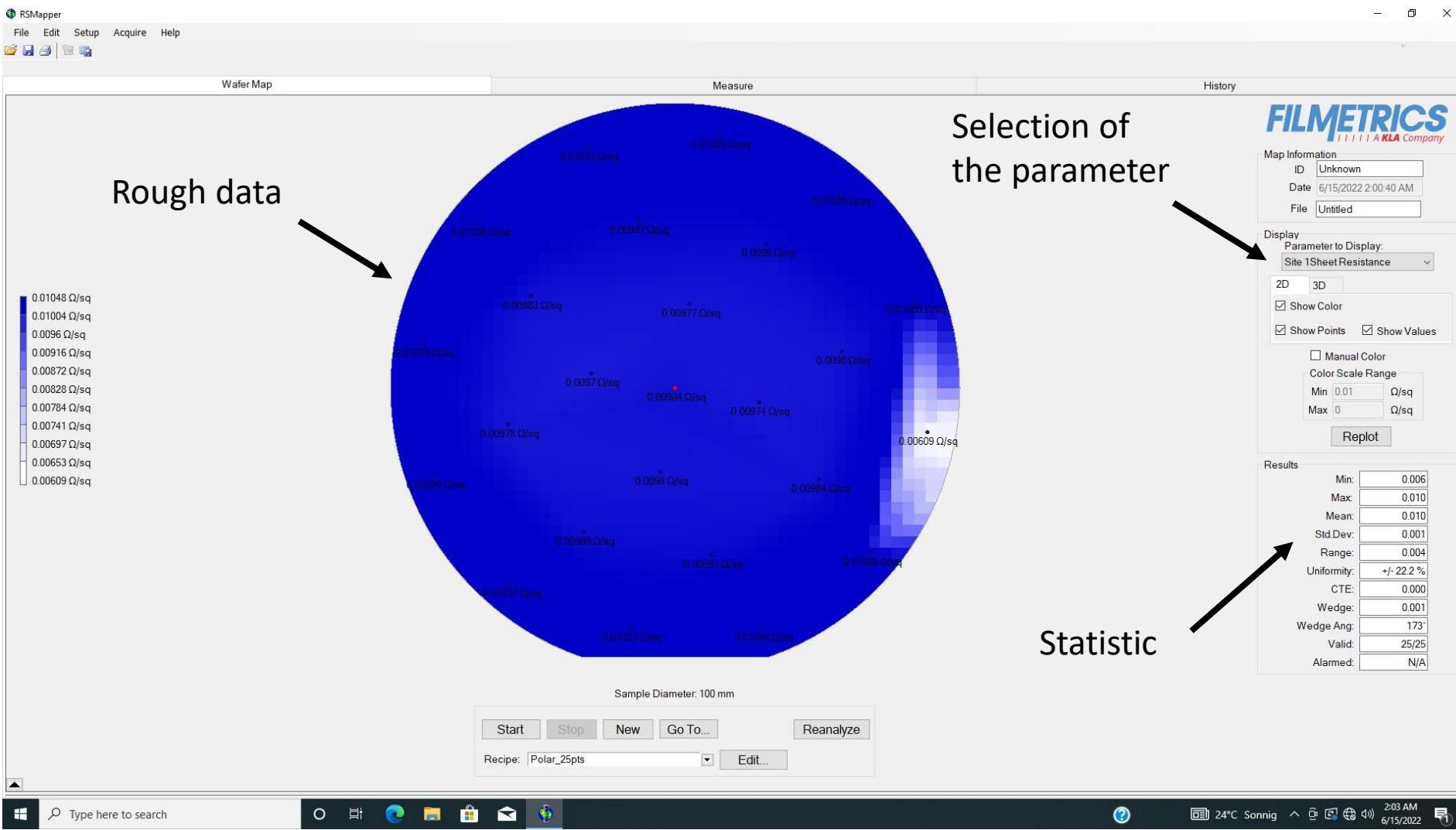

Visualization of the measurements on the “History” tab

Statistic

Rough data

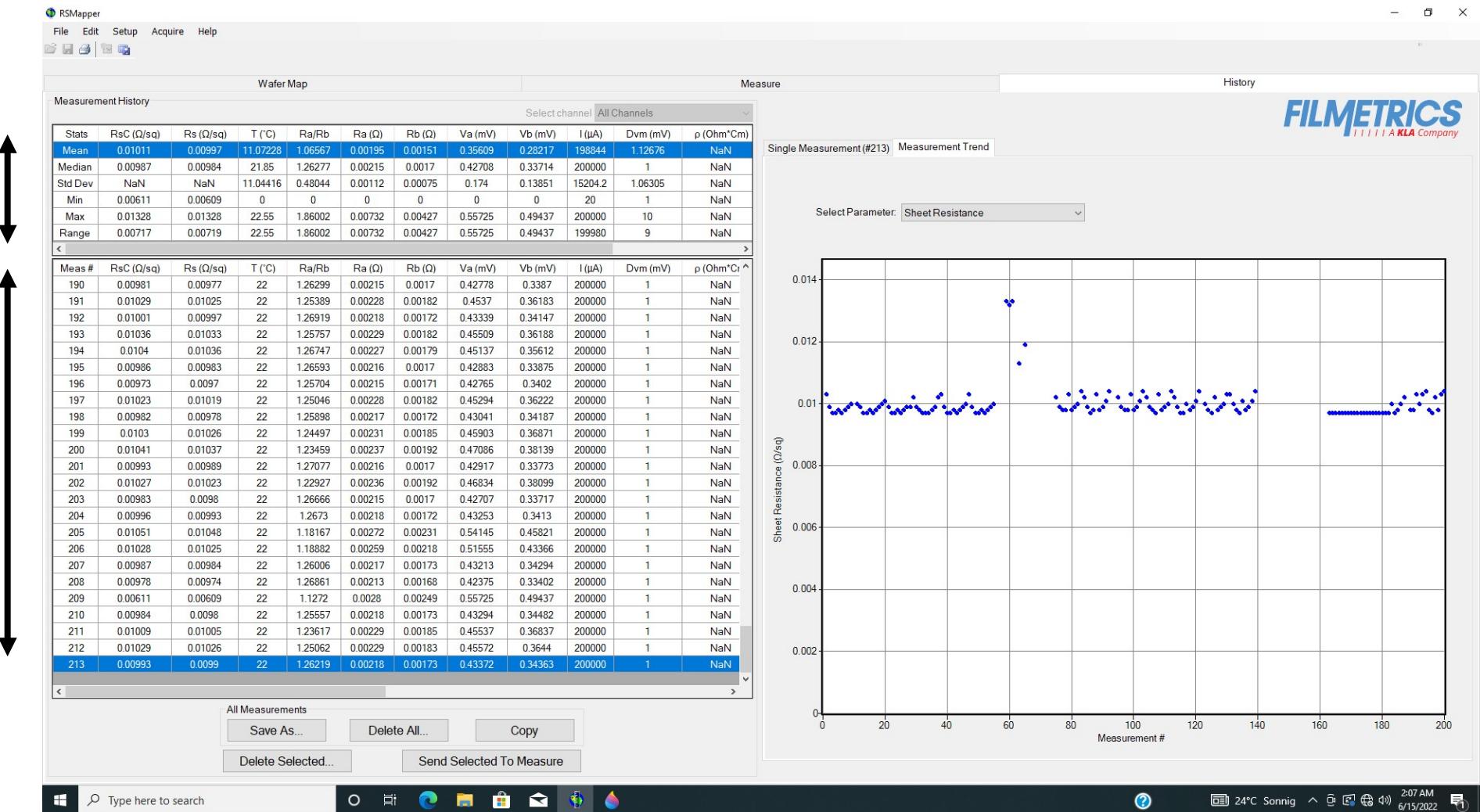

11. Once the measurement is done, click the «Go To» button




11

12. Click “Go to Unload”

13. Take your wafer / sample

14. Close the software

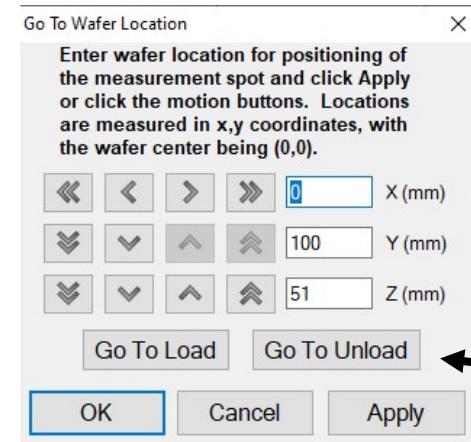

12