



Alliance Concept EVA760 Uniformity Masks Installation

1. 450 mm

• After installing the dome, pull the mask upwards until it is clipped together horizontally.





Figure 1: Left : Mask 450 mm in down position. Right : Mask 450 mm in up position, ready for the evaporation.





2. 350 mm and 250 mm

• First, without the dome installed, put the 450 mm Mask in the horizontal position (see point 1) and remove the pin. Be careful when removing the pin to not damage / bend it.





 Put the 450 mm mask back to its down position, rotate it fully to the left and place the pin back.







 Untight and remove the two the screws, place either the 250mm or the 350 mm and tight the 2 screws fully back in. The 250 mm mask goes on the bottom slot, the 350 mm one on the top slot. The mask is now ready.



Figure 2: Left : 250 mm Mask screwed in place. Right : 350 mm Mask screwed in place.

• After the evaporation, follow the protocol in the inverse order to dismount the mask. Please tight the screws fully back in to avoid any material to deposit on the thread.

