

Sort your BIO waste (P1 only) STI-IN 2023

Security – Stability

Temperature, pH, swelling, evaporation, smell, container condition

Liquid
Waste

Solid
Waste

Biological P1
Inactivated with
Bleach or Virkon
18 01 06

Needles
sharp objects

Non-genetically
modified and non-
infectious
organisms -
gloves,
contaminated
papers, tips,
empty tubes and
flasks, fixed
tissues, etc.

Material
contaminated with
GMOs AND
possibility to
dispose of sharp
objects (pipette
tips, serological
pipettes, tubes,
etc.; EXCEPT
needles and
scalpel blades)



Let the bio solution et
virkon or bleach degas for
at least **1 day** in a chemical
hood with the cap slightly
open to release the gas.

Needles
disposal
container
18 01 01

Biological waste
to be deposited in a yellow
container or in a designated room.

Biological waste to be deposited in an STI
collection center

